





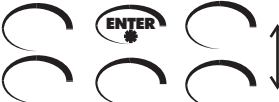












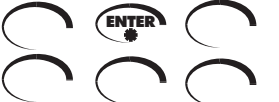


Quick Start Guide

1.	Press  to turn on Spectro. MAIN MENU will appear.		<pre> MAIN MENU 12:45 * CALIBRATE WL PROGRAMMED TESTS %T/ABS </pre>
2.	Press  or  to scroll until the * is next to PROGRAMMED TESTS.		<pre> MAIN MENU 12:45 CALIBRATE WL * PROGRAMMED TESTS %T/ABS </pre>
3.	Press  to select PROGRAMMED TESTS. PROGRAMMED TESTS menu will appear.		<pre> PROGRAMMED TESTS * SEQUENCE 1 SEQUENCE 2 SEQUENCE 3 </pre>
4.	Scroll to ALL TESTS with  button.		<pre> PROGRAMMED TESTS SEQUENCE 2 SEQUENCE 3 * ALL TESTS </pre>
5.	Press  to select ALL TESTS. ALL TESTS menu will appear.		<pre> ALL TESTS * 1 Aluminum 2 Alkalinity-UDV 3 Ammonia-N L F </pre>
6.	Press  or  to scroll to desired test.		<pre> ALL TESTS 76 pH TB 77 Phenol * 78 Phosphate L </pre>
7.	Press  to select test. Menu for desired test will appear.		<pre> 78 Phosphate L SCAN BLANK * SCAN SAMPLE END 880NM </pre>
8.	Insert blank, press  to scan blank.		<pre> 78 Phosphate L SCAN BLANK * SCAN SAMPLE END 880NM </pre>
9.	Insert reacted sample. Press  to scan sample. Result will appear.		<pre> 78 Phosphate L 74.3%T 0.1297A 0.27PPM PRINT PRESS ENTER </pre>